AFI present
THE GENERATION

KATRINA

BLOOD BANK & HIGH CAPACITY CENTRIFUGE

From Experience comes Excellence

2022

UNIQUE FRENCH MANUFACTURER

AFI Centrifuge, the only French manufacturer based in Château-Gontier in FRANCE, has been designing and manufacturing innovative centrifuges since 2009. AFI centrifuges have been designed to meet the needs and requirements of medical analysis laboratories, research laboratories, veterinary laboratories, hospitals, etc.

Offering a wide variety of rotors and accessories, the user will have at his disposal all the solutions to meet the required protocols. Available in both ventilated and refrigerated versions, AFI centrifuges are designed to be used with all types of tubes, vials and bottles available on the market.

With more than 20 years of experience in the field of biomedical engineering, electronics and electromechanics, AFI presents today a wide range of versatile and innovative centrifuges

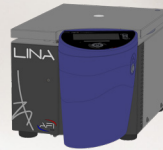
Philippe LE GUYADER

A stylized signature of Philippe Le Guyader in blue ink.

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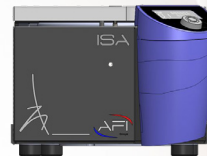
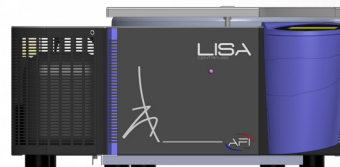
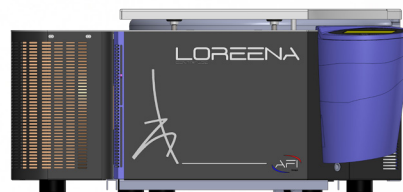
REGINA
High Capacity Centrifuge
12 L

LIA

Rotor 24 micro tubes
15 000 Rpm/ 21 380 x g
Ventilated & Refrigerated
8 rotors

LINA

Rotor 48 micro tubes
Rotor 6x 50mL
Rotor 6 Micro plates
23 000 Rpm/ 39 625x g
Ventilated & Refrigerated
11 rotors

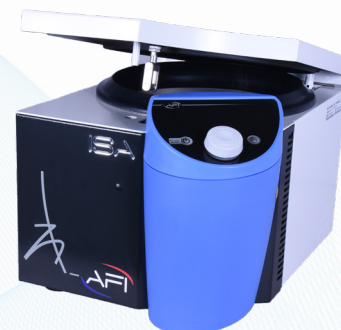
SIRENA

24-36 tubes 5/7 mL
6x 100mL
18 000 Rpm/ 31514 x g
Refrigerated
Interactive display
Touch screen, 99 programs,
17 rotors

ISA

40 tubes 5/7 mL
6700 Rpm/ 6324 xg
Ventilated & Refrigerated
Touch Screen, 99 programs,
11 rotors

Micro
Research centers / Veterinaire

Performance
Research centers/ Universities

Compact
Small laboratories

High Capacity
Blood bank / Industry

KATRINA

12x Bags 500mL / 4x 2 Liters

4900 Rpm/ 7301 x g

Refrigerated Floor

Interactive communication, Touch screen,
Integral, 6 Rotors

LOREENA

108- 200 tubes 5/7 mL / 12x 50 mL

16000 Rpm/ 28 335x g

Ventilated & Refrigerated & Floor

Interactive communication,
Click Spin, Touch Screen,
15 Rotors

LISA

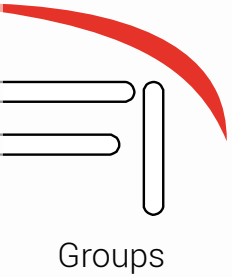
60-92 tubes 5-7 mL/ 8x100mL

18 000 Rpm/ 31514 x g

Ventilated & Refrigerated

Interactive Communication, Click Spin, Touch
Screen, 18 Rotors

Routine
Hospitals/ Medical analysis Laboratories

KATRINA

BLOOD BANK Centrifuge

NEW GENERATION

The new generation of centrifuge for your blood bank center KATRINA benefits from the experience of AFI, it will easily and safely replace your old centrifuges. Katrina meets all your validated protocols for separating plasma, Buffy coat, red blood cells.

QUALITY THAT SPEAK FOR ITSELF

The «IDEM» application coupled with PFC (Power Factor Control) technology ensures a reproducibility of your cycles making Katrina an ideal centrifuge for simplified blood processing productivity

INTERACTIVE COMMUNICATION

With its large «Smartphone» touch screen colour display, KATRINA ensures interactive communication between you and the machine. Backlit display for better visibility Programming is greatly facilitated by direct access to all cycle parameters in a single window. Compatible with gloved hands.


SECURITY

User access as password protection of the protocols ensures increased security and efficiency

Centrifuge	Description
AFI-C600R-2P-B12	KATRINA Blood Banking Centrifuge 12x450/500 mL
AFI-C600R-2P-B6	KATRINA Blood Banking Centrifuge 6x550/600 mL
Specifications	Swing-out rotor
Max Speed	4900rpm
Max RCF	7301 g
Max Loading	6x1000ml
T° Range	-20°C/+40°C
Dim. (HxWxD)	95x81x96cm
Pack dim.	114x89x109cm
Weight	352kg

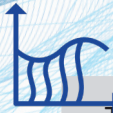The high performance centrifuge

IDEM APPLICATION

The IDEM application guarantees a perfect reproduction of the cycle. Regardless of the environment (rotor change, partial bag loading, tension, machine age) the IDEM application guarantees identical results for all your cycles.

INTERNATIONAL AUTOMATIC ELECTRICAL CORRECTION

International automatic correction of the electrical network. Guarantees stability of the machine's performance.

INTERACTIVE DISPLAY

Its large touch screen, ensures simple communication between user and device.
Backlight display for better visibility
One screen that grants direct access to all cycle parameters, also compatible with gloves

Its major assets

6 reasons

to choose it

Multi-user, password-protected access control for protocol protection

SECURITY

Centrifuge's electronic functions integrated into a front central unit. Plug & play system guaranteeing a fast and efficient intervention

THE BLUE BOX

You can set up your own programs, name it, lock it, choose slope rate, set up your own user preferences.

CUSTOMIZABLE PROGRAM

KATRINA

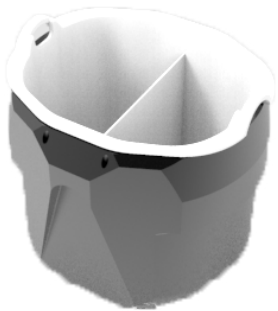
KATRINA, AFI's Topnotch Centrifuge

Designed to meet all needs related to blood treatment and optimized to reach the highest level of performance, while remaining ergonomic and user-friendly.

KATRINA, the New Generation in Centrifuge for your Transfusion Centers

Optimally designed by a group of French experts, KATRINA is a safe and easy alternative to your old centrifuges. Katrina meets validated protocol and international standards.

KATRINA, Quality that Speaks for Itself

The « IDEM » application combined with PFC (Power Factor Control) technology ensures cycle reproducibility, making KATRINA the ideal centrifuge for simple and productive blood treatment.

NEW GENERATION

KATRINA, AFI's Savoir-Faire

Modelled after AFI's latest technical practices, KATRINA truly represents state-of-the-art French design.

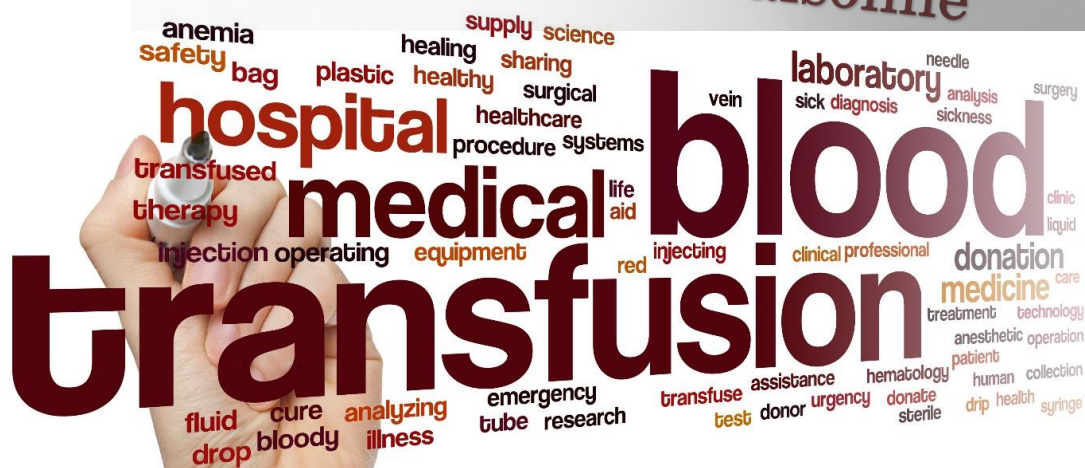
- AFI designs and manufactures its own motors
- AFI designs and manufactures its own circuit boards
- AFI designs and manufactures its own cooling units.
- AFI designs and manufactures its own parts
- Its innovative suspension and noise reduction system make the AFI centrifuge the most reliable of its kind.

All AFI centrifuges are manufactured in Mayenne (France), a longtime leader in French centrifugation, and sealed with the "Made in France » quality stamp of approval


KATRINA

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
AFI RX6-1000-12

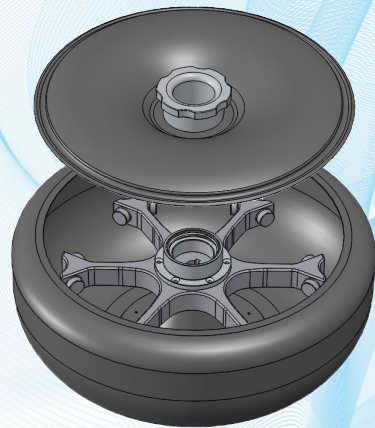Blood Bags Adapters selection guide

Model	R
Max Capacity	6x 1L
Angle	90°
RPM Max	4 300
ACR Max	5 767

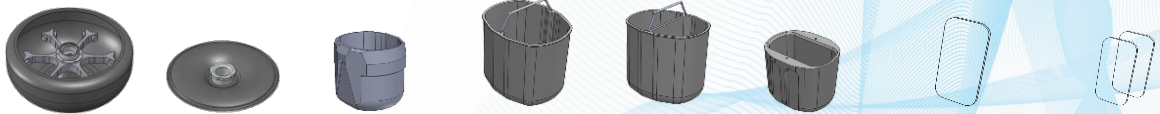Accessories

Ref.	AFI-RX6-1000-12	AFI-LRX6**	AFI-BO6-1000*	AFI-LI6-1000*	AFI-LIS6-1000*	AFI-6BD1000	AFI-6TBB	AFI-WLI750*
		* Set of 6		** Recommended for use above 3800 rpm				* Set of parts from 10 to

Blood Bag	450 mL	500 mL	550 mL	600 mL	750 mL
Type	Double to Quadruple	Quintuple Quintuple + F	Quadruple Quintuple	Simple Double	Simple
Bags per adapter / rotor	1 6	1 6	1 6	1 6	1 6

Adapter

Hole LWH (mm)	1 Cavity Liner 143x116x143				
Radius	279				
Max charge (gr)	1800				
Ref.	AFI-LI6-1000				

Blood Bags	350-450 mL	450 mL	500 mL	1L
Type	Simple Double	Quadruple	Triple	Simple
Bags per adapter / rotor	2 12	2 12	2 12	1 6


Adapter

Hole LWH (mm)	2 Cavities Liner 69x116x143			70 x 130 x 120
Radius	279			270
Max charge (gr)	1800			
Ref.	AFI-LIS6-1000			AFI-6BD1000


AFI RX6-1000-6

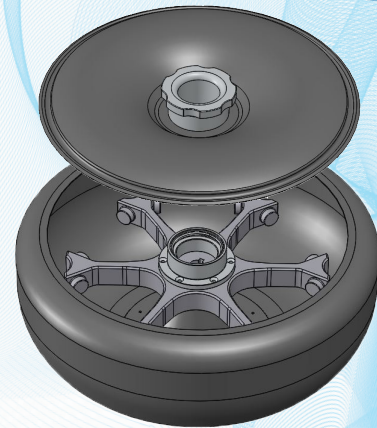Blood Bags Adapters selection guide

Model	R
Max Capacity	6 x 950ml
Angle	90°
RPM Max	4 900
ACR Max	7 301

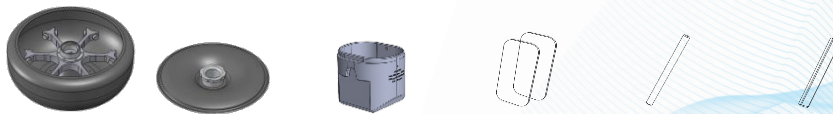Accessories


Ref.	AFI-RX6-1000-6	AFI-LRX6*	AFI-BO6-750	AFI-WLI750*	AFI-STICK7	AFI-STICK14
------	----------------	-----------	-------------	-------------	------------	-------------

** Recommended for use above 3800 rpm

* Set of parts from 10 to 40gr

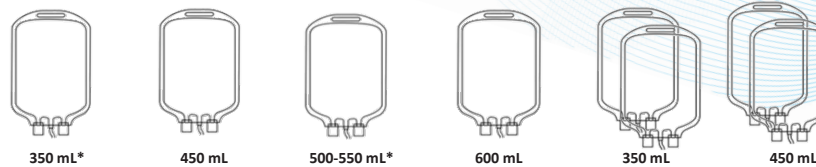Blood Bags

Type	Double / Triple Quadruple + Filtre	Double / Triple Quadruple + Filtre	Double / Triple Quadruple + Filtre	Simple Buffy coat run
Bags per adapter / rotor	1 6	1 6	1 6	2 12

Adapter

Hole LWH (mm)				
Radius	272	269	269	269
Max Charge (gr)	1500			
Ref.	AFI-LI6-750	AFI-LI6-750-600	AFI-LI6-750-600	AFI-LI6-750-900

* We recommend the use of a silicone spacer to wedge the pocket

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KATRINA

Centrifuge HIGH PERFORMANCE

The **KATRINA** High Performance has been designed to meet the bioprocessing methods.

KATRINA High Performance is available with a choice of 2 rotors : 6x1L or 4x2L to meet all your needs for high throughput sample processing.

PROVEN QUALITY

The «IDEM» application coupled with PFC (Power Factor Control) technology ensures reproducible cycles, making **KATRINA** an ideal centrifuge for simplified productivity.

INTERACTIVE COMMUNICATION

With its large «Smartphone» touch screen colour display, **KATRINA** ensures interactive communication between you and the machine. Backlit display for better visibility. Programming is greatly facilitated by direct access to all cycle parameters in a single window. Compatible with gloved hands.

SECURITY


User access as password protection of the protocols ensures increased security and efficiency

Centrifuge	Description
AFI-C600RE-2P-H6	KATRINA Haute Performance centrifuge
AFI-C600RE-2P-H8	KATRINA Haute Performance centrifuge
Specifications	Swing-out rotor
Max Speed	4300 / 4400 rpm
Max RCF	5809 / 5931g
Max Loading	6x1L / 4x2L
T° Range	-20°C/+40°C
Dim. (HxWxD)	95x81x96cm
Pack dim.	114x89x109cm
Weight	352kg

Swinging bucket rotors

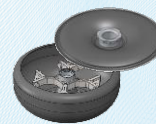Accessories selection guide

KATRINA ACCESSORIES	AFI-RX6-1000-H	AFI-RX6-1000-D	AFI-RX4-2000
---------------------	----------------	----------------	--------------

KATRINA ACCESSORIES	AFI-BR6-1000	AFI-BXDW6	AFI-BO6D-1000*	AFI-BR4-2000
Max Capacity (ml)	6x 1L	48 Microplates	6x 1L	4x 2L
Max Charge (gr)	1400	1000	1800	2400
Max Speed (rpm)	4 300	4 300	3 600	4 400
Max RCF (g)	5 747	5 519	4 289	5 931
Cavity dim (ØxD) or (WxLxD)	Ø102x135	90x132x114	117x144x152	Ø124x144
Lid	✓	✓		


Bottles	ml	Ø x L				
	150 FALCON	61x118	6	-	-	4
	200 FALCON	61x138	6	-	-	4
	200 AFI-B-PPCO200C	56x118	6	-	-	4
	200 NUNC/NALGENE	60x130	6	-	-	4
	200 AFI-B-PPCO200F	56x112	6	-	-	4
	250 CORNING	60x163	6	-	6*	4
	250 AFI-B-PPCO250F	62x122	6	-	6*	4
	290 AFI-B-PPCO290F	62x137	6	-	6*	4
	400	81x136 84x134	6	-	6*	4
	500 CORNING	96x148	6	-	6*	4
	500 AFI-B-PPCO500F	98x114	6	-	6*	4
	625 AFI-B-PPCO625F	92x130	6	-	6*	4
	750 AFI-B-PPCO750F	98x152	6	-	6*	4
	750 AFI-B-PPCO750FB	100x130	6	-	6*	4
	950 AFI-B-PPCO950FB	100x154	6	-	6*	4
	1L AFI-B-PPCO1000FB	100x168	6	-	6*	4
	1L AFI-B-PPCO1000F	98x175	6	-	6*	4
	1L NALGENE	94x217 94x197	-	-	6*	4*
	2L AFI-B-PPCO2000F	125x214	-	-	-	4
	Micro Plates		-	48	-	-
	Deep Well		-	12	-	-

*UNSEALED


AFI RX6-1000-H

Adapters selection guide

Model	R
Max Capacity	6x 1L
Angle	90°
RPM Max	4 300
ACR Max	5 499

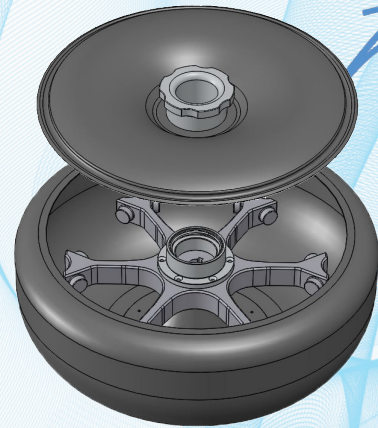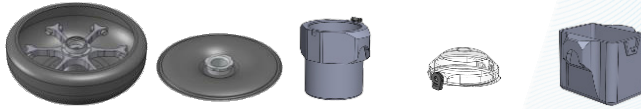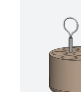
Accessories

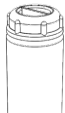
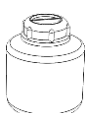
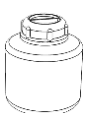
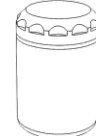
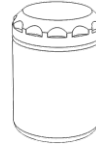
Ref.	AFI-RX6-1000-H	AFI-LRX6**	AFI-BR6-1000*	AFI-LR6-1000	AFI-BXDW6*
------	----------------	------------	---------------	--------------	------------

* Set of 6
** Recommended for use above 3800 rpm

Tubes									
Capacity (ml)	15	15	30	30	30	50	50	50	50
∅ x L (mm)	17x120	17x120	25x110	25x110	25x95	29x115	29x115	29x115	29x107
Tubes per adapter / rotor	13 78	13 78	7 42	7 42	8 48	5 30	5 30	5 30	7 42
Adapter									
Hole ∅ x L(mm)	18x80	18x80	26x60	26x60	26,5x70	30x70	30x70	30x70	30x70
Radius	278	278	271	278	276	271	271	278	276
Max Charge (gr)	1400	1400	1400	1400	1400	1400	1400	1400	1400
Ref.	AFI-6T15C-RB	AFI-6T15ST-RB	AFI-6T30S-RB	AFI-6T30C-RB	AFI-6T30AOR-RB	AFI-6T50S-RB	AFI-6T50CS-RB*	AFI-6T50C-RB	AFI-6T50AOR-RB

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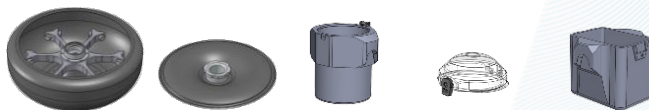
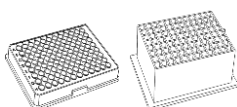
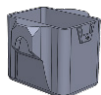
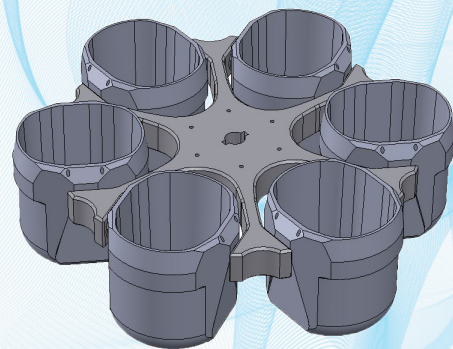
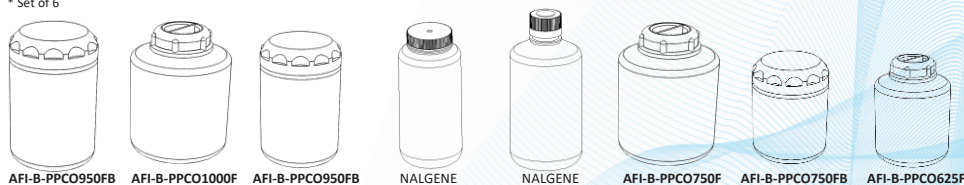
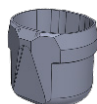
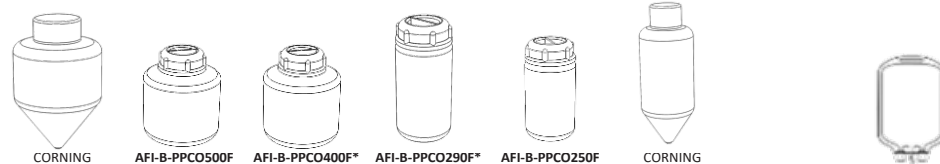
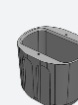
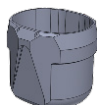
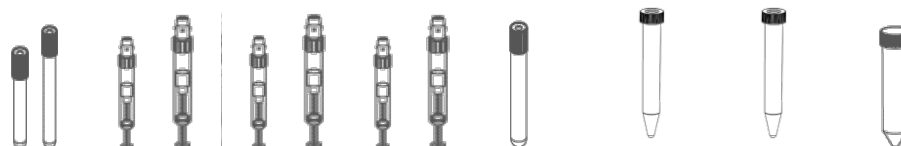
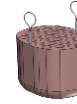
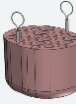
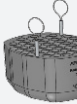
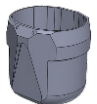
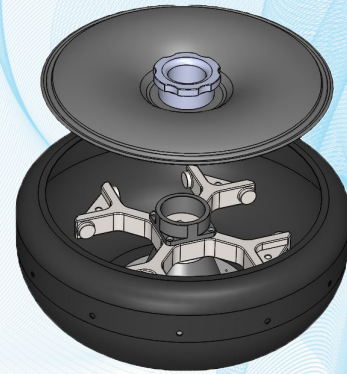
KATRINA

Centrifuge AUTOMATION

The centrifuge **KATRINA** Automation have been designed to meet the automation systems used in laboratories. **KATRINA** Automation is compatible with a wide range of racks that can be centrifuged directly in dedicated buckets. An undeniable time saving in the analysis.

PROVEN QUALITY

The «IDEM» application coupled with PFC (Power Factor Control) technology ensures reproducible cycles, making **KATRINA** an ideal centrifuge for simplified productivity.

INTERACTIVE COMMUNICATION

With its large «Smartphone» touch screen colour display, **KATRINA** ensures interactive communication between you and the machine. Backlit display for better visibility. Programming is greatly facilitated by direct access to all cycle parameters in a single window. Compatible with gloved hands.

SECURITY

User access as password protection of the protocols ensures increased security and efficiency.


Centrifuge	Description
AFI-C600RE-2P-AUTO	KATRINA Automation
Specifications	Swing-out rotor
Max Speed	3300 rpm
Max RCF	2313g
Max Loading	3,8kg
T° Range	-20°C/+40°C
Dim. (HxWxD)	95x81x96cm
Pack dim.	114x89x109cm
Weight	352kg

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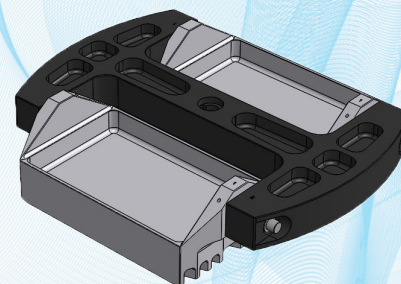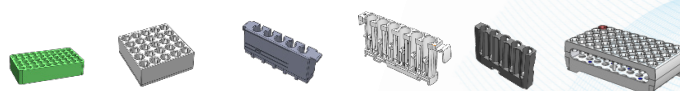
KATRINA

Centrifugeuse VETERINARY

The **KATRINA** Veterinary centrifuge has been designed to meet the intense analysis periods of veterinary laboratories.

Its dedicated accessory and its automatic tube uncapping system during centrifugation offer an undeniable time saving to laboratories during the prophylaxis period for example.

192 tubes can be centrifuged and uncorked at the same time.

PROVEN QUALITY

The «IDEM» application coupled with PFC (Power Factor Control) technology ensures reproducible cycles, making **KATRINA** an ideal centrifuge for simplified productivity.

INTERACTIVE COMMUNICATION

With its large «Smartphone» touch screen colour display, **KATRINA** ensures interactive communication between you and the machine. Backlit display for better visibility. Programming is greatly facilitated by direct access to all cycle parameters in a single window. Compatible with gloved hands.

SECURITY

User access as password protection of the protocols ensures increased security and efficiency.


Centrifuge	Description
AFI-C600RE-2P-V	KATRINA Veterinary centrifuge
Specifications	Swing-out rotor
Max Speed	3300 rpm
Max RCF	2252g
Max Loading	192 tubes
T° Range	-20°C/+40°C
Dim. (HxWxD)	95x81x96cm
Pack dim.	114x89x109cm
Weight	352kg

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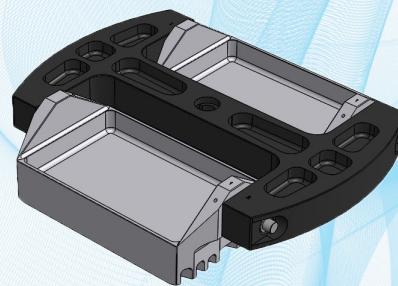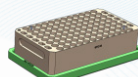
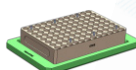
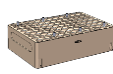
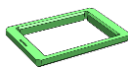
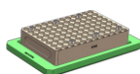
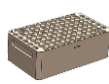
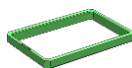
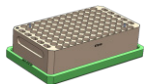
e-mail: contact@afigroups.com

*Scan me to discover
AFI catalogs*

* If necessary or useful without affecting the essential characteristics of a model, the technical and stylistic characteristics may be modified without prior notice from the manufacturer.